Data Capture revision history

Version 6.0.1 4/1/2021

Added function to update the time/date chip on all programmed units. This is great for updating for Day Light Savings.

Version 6.0.1 3/10/2021

The big change is writing sensor data to it's own individual database ie. tblSensorDatalD_xxx.fic. This was done to speed up reports.

Added ability to disable email reminder

Version 5.4.10

Added query return all sensor data for each unit to add these reading to the daily status email.

Added ability to add contact name for each phone number. This is a feature requested by a customer.

Version 5.4.10 1/30/2020

Added support for M408/M409 CO2 monitor.

Version 5.4.10 11/3/2020

Corrected code for downloading log for model VM608E.

Version 5.4.10 10/21/2020

Added code to show IP address in Temperature Guard Units tree.



Version 5.4.9 9/22/2020

Corrected the units displayed in the Hi-Lo report for humidity sensor. Increased font size on all menus.

Version 5.4.8 4/28/20

Added button to check and repair and re-index all database files.

Version 5.4.8 1/2020


Added button to delete unsent alarm emails to the configuration page.

Moved Save, log file buttons to menu bar.


Data Capture revision history

Rev 5.4.7 Compiled and RELEASED 11/8/2019

A local hospital customer requested the ability to edit all unit email addresses at once. Added menu item under Edit

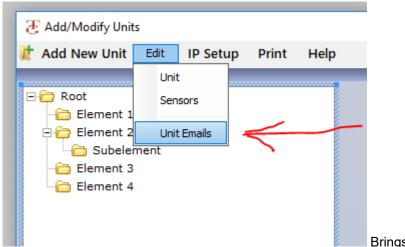

Brings up below screen

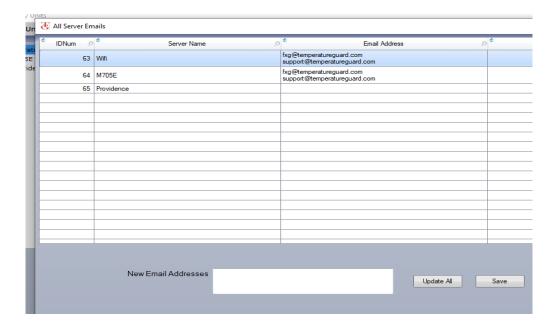

Rev 5.4.6 10/28/2019

Customer requested the Alarm Email Reminder options include 5 minutes. The available reminder times are changed depending on the update rate.